Aseptic Sampling Valves

Sanitary
flow
equipment

Tipo **VPAK**

Pres campion asettica

CARATTERISTICHE

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
STERILE SAMPLING

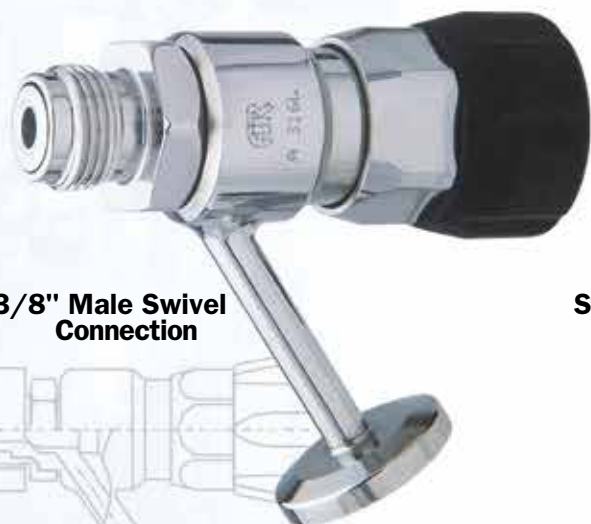Developed for taking samples in a sterile way, the special configuration assures effective cleaning (CIP) and sterilization (with saturated steam) of product contact surfaces of valve and seating plug with the valve in closed position. Once installed, the sealing plug shuts off flush with the inlet port connection with a zero dead leg.

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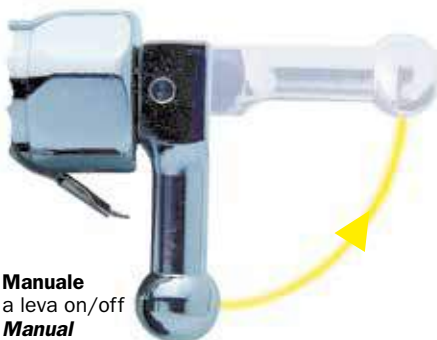
Manuale
a leva on/off
Manual
stainless steel on-off lever

Pneumatico inox on/off
Air operated stainless steel air-spring actuators

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DOCUMENTATION

All valves are fully traceable for validation processes. Each valve is accompanied by material test certificates EN 10 204 3.1B 3A standards compliance certificates, FDA regulations compliance certificates, and tapes detailing the valve interior surfaces Ra.

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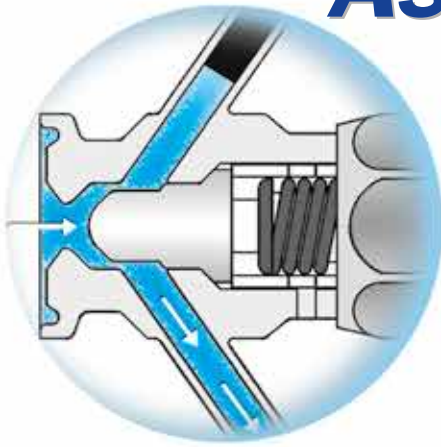
IMPORTANT!

For reliable performance there must be a safety margin between the spring load of the seating plug and the maximum pressure in the product line.

Higher pressures will force open the valve.

Please indicate the maximum pressure of your applications.

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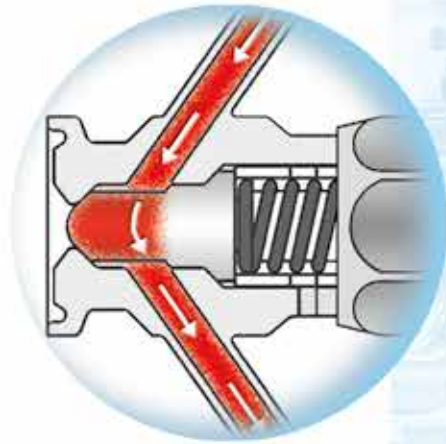
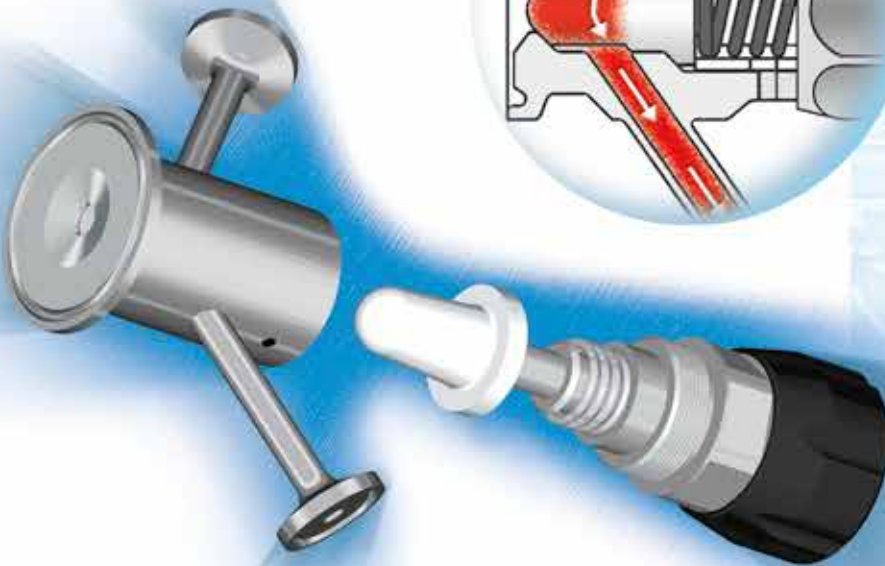
Valvola chiusa (fase di CIP/steriliz.)
Valve closed (CIP/steril. phase)

VPA-VPK Ordering information

To specify the part completely, start with the product description and select the additional options as shown below :

Valve	Option	Inlet hole	Inlet	Sampling	CIP/SIP	Diaphragm	Actuator	Sensor	finish	material
VPK	-	05	AG	PB	AG	T	MC		-11	A

Valve	Inlet Hole
VPK	05 5 mm
VPA	07/10/13 7 / 9,4 / 12,7 mm
VPZ	05/07/10/13
on request	(*) 7mm Silicone only

Option
- Standard Safe Supply
A ATEX execution

Inlet Connection	VPK	VPA
clamp	1/2" AG	•
	3/4" AH	•
	1" AJ	•
	1 1/2" AM	•
	2" AN	•
butt weld	Ø19mm QC	•
	Ø30mm QF	•
BSP screw	1/4" LC	•
	3/8" LE	•
	1/2" LG	•
Ingold	Ø25 ZA	•
	flanged S.	•
on request		

Sample - CIP/SIP	VPK	VPA
clamp	1/2" AG	•
	3/4" AH	•
to be weld	Ø6,25x1,24mm RB	•
	Ø12,7x1,65mm BG	•
hose barb	Ø6,25mm PB	•
	Ø12,7mm PF	•
quick connect	ZG	•
on request		
Sensor		
11	Inductive PNP 12/24VdC	
on request		

Diaphragm	
T TFM PTFE	
X Silicone	
Actuator	
MC	Manual wheel
ML	Manual lever on/off
MR	Lever auto return
PN	Pneumatic s.e./n.c.
PL	Pneumatic + lever

Surface Finish VPK	
11 ≤ 0,5 µm	
51 ≤ 0,5 µm + Electropolish	
Surface Finish VPA	
71 ≤ 0,5 µm	
41 ≤ 0,4 µm + Electropolish	
on request	
Materials	
A	AISI 316L-1.4404
B	1.4435
C	C22 Hastelloy
D	C276 Hastelloy
on request	

Aseptic sampling valve

Type **VPA**

CHARACTERISTICS

The **AR** sampling valve range has been designed for sterile sampling applications in the Chemical and Pharmaceutical industries. Its compact design allows easy installation on small vessels or process lines.

FULLY AUTOCLAVABLE,

Valve body and head are made out of certified 316L bar stock.

Damage to the seating plug and unexpeted material in the outlet are indications of overpressure. A leak detector hole is placed immediately behind the outlet.

On request the valve can be equipped with a CIP system.

SURFACE FINISH

External mechanical "mirror" polish
Internal mechanical polish Ra < 0.4 µm

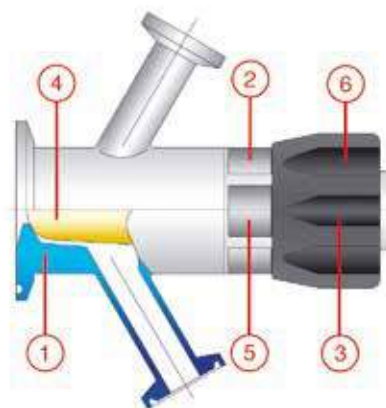
VALVE BODY connections available in:

- weld ends
- clamp ends

Other types available upon request.

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